

# ECLIPSE Ti2

Inverted Research Microscope



Shedding New Light On **MICROSCOPY** 

# See More Than Before

## Leading platform for advanced imaging

The ECLIPSE Ti2 delivers an unparalleled 25mm field of view (FOV) that revolutionizes the way you see. With this incredible FOV, the Ti2 maximizes the sensor area of large-format CMOS cameras without making compromises, and significantly improves data throughput.

The Ti2's exceptionally stable, drift-free platform is designed to meet the demands of super-resolution imaging while its unique hardware-triggering capabilities enhance even the most challenging, high-speed imaging applications. Furthermore, the Ti2's unique, intelligent functions guide users through imaging workflows by gathering data from internal sensors, eliminating the possibility of user errors. In addition, the status of each sensor is automatically recorded during acquisition, providing quality control for imaging experiments and enhancing data reproducibility.

In combination with Nikon's powerful acquisition and analysis software, NIS-Elements, the Ti2 is a total innovation in imaging.


## **Contents**

Groundbreaking FOV	P4	Intuitive operation —	P12
High-quality Nikon optics	P6	System expandability	P14
Focus perfected —	P8	NIS-Elements	P16
Assist Guide —	P10	System diagrams	P18



Motorized and intelligent model for advanced imaging applications. Compatible with PFS, auto correction collar, and external phase contrast system. The base of choice for live-cell imaging, high-content applications, confocal and super-resolution.


Manual model with imaging capability for laser applications. Intelligent features provide interactive guidance through imaging workflows and automatically detect microscope status.

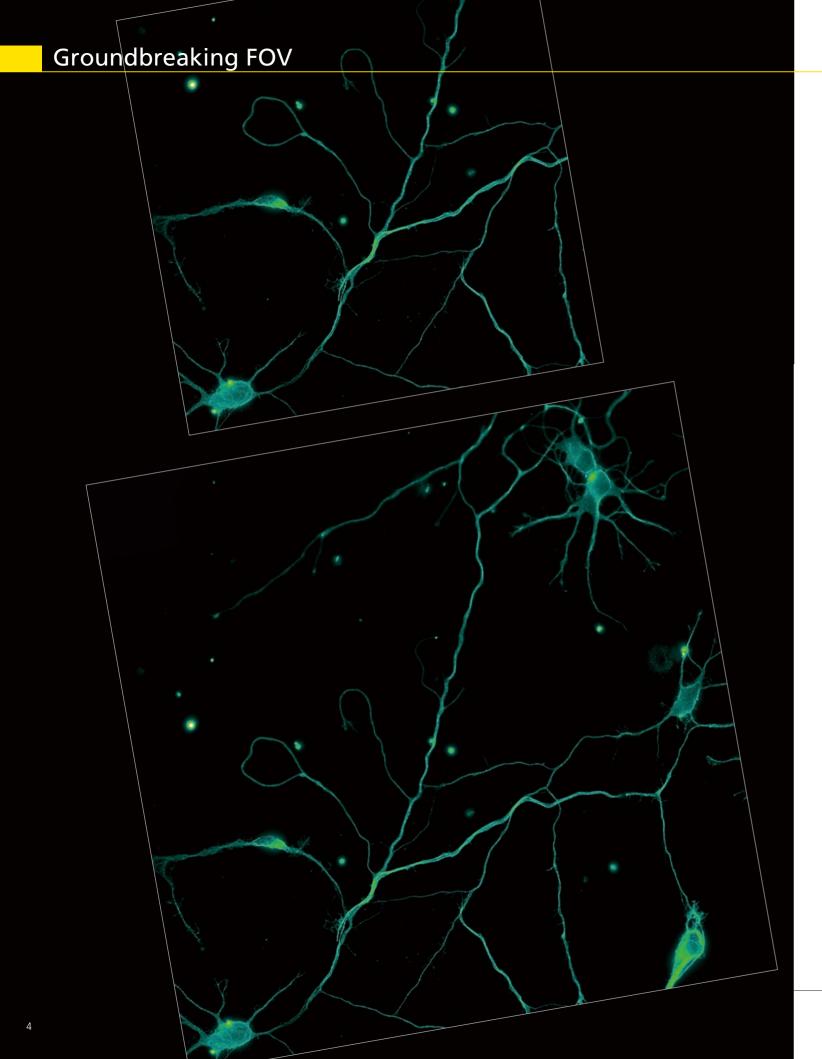


As research trends evolve towards large-scale, systems-level approaches, there is an increasing demand for faster data acquisition and higher throughput capabilities. Development of large-format camera sensors and improvements in the data processing capabilities of PCs have facilitated such research trends. The Ti2, with its unprecedented 25mm field of view, provides the next level of scalability, enabling researchers to truly maximize the utility of large-format detectors and future-proof their core imaging platform as camera technologies continue to develop at a rapid pace.

### Bright illumination over a wide area

High-power LEDs deliver bright illumination across the Ti2's large field of view, ensuring clear, consistent results from demanding applications such as high-magnification DIC. Incorporation of a flyeye lens design provides uniform illumination from edge to edge for quantitative high-speed imaging and seamless tiling of images in stitching applications.

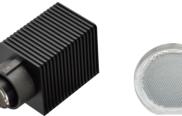

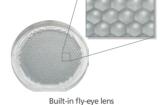

A compact epi-fluorescence module designed for large FOV imaging provides high transmittance across a broad spectrum, including UV. Large diameter fluorescence filters with hard coatings deliver large FOV images with a high signal-to-noise ratio.



Eni-fluorescence module Large diameter fluorescence filter cubes

## Large diameter observation optics

The diameter of the observation light path has been enlarged in order to achieve a field number of 25 at the imaging port. The resulting large FOV is capable of capturing approximately double the area of conventional optics, enabling users to gain maximum performance from large-format sensors such as CMOS detectors.



High-power LED illuminator





Imaging port with large 25 field number

## **Objectives for large FOV imaging**

Objectives with superior image flatness ensure high quality images from edge to edge. Utilizing the maximum potential of the OFN25 objective significantly accelerates data collection.



## Cameras for large-volume data acquisition

Nikon's FX-format F-mount cameras Digital Sight 10 and Digital Sight 50M are equipped with CMOS image sensors optimized for research use, originally developed for professional D-SLR cameras. This allows high-speed and high-sensitivity live-cell imaging, enabling the best use to be made of the



D-SLR camera technology optimized for microscopy








Digital Sight 50M

Ti2's large FOV.

Neuron culture stained for microtubules (Alexa Fluor® 488), captured with CFI Plan Apochromat Lambda 60XC objective and DS-Qi2 camera Conventional FOV on top and new Ti2 FOV on bottom

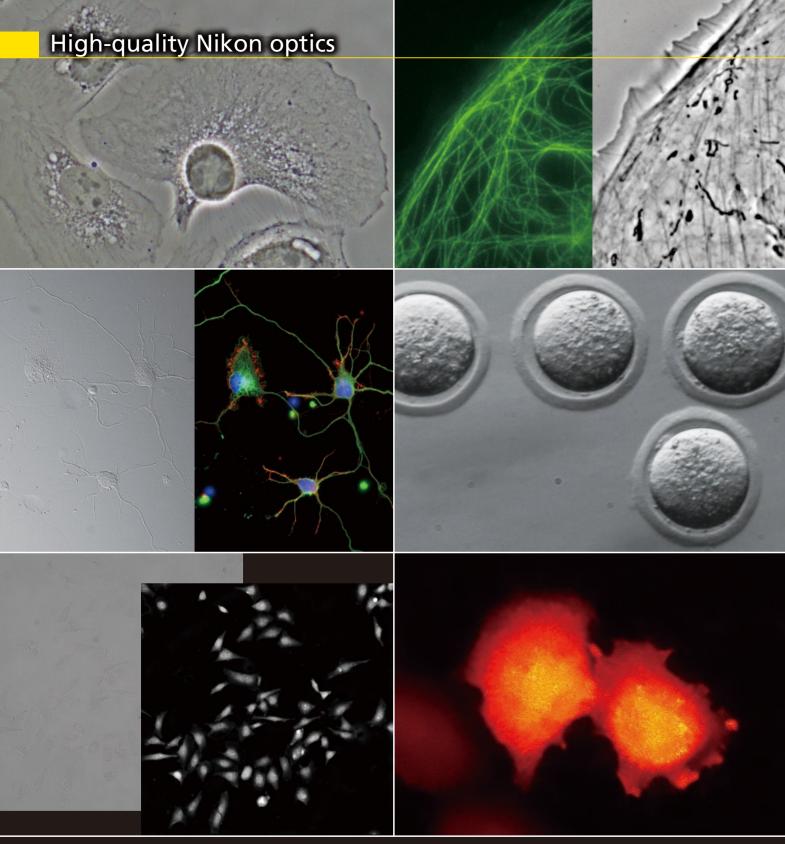

Apodized phase contrast image: BSC-1 cells captured with CFI S Plan Fluor ELWD ADM 40XC objective

DIC and epi-fluorescence images: 25mm FOV image of neurons (DAPI, Alexa Fluor® 488, Rhodamine-Phalloidin), captured with CFI Plan Apochromat Lambda 60XC objective and DS-Qi2 camera Photo courtesy of Josh Rappoport, Nikon Imaging Center, Northwestern Univ.; Sample courtesy of S. Kemal, B. Wang, and R. Vassar, Northwestern Univ.

Brightfield and Volume Contrast images: HeLa cells captured with CFI S Plan Fluor ELWD 20XC objective

Epi-fluorescence and external phase contrast images: PTK-1 cells labeled with GFP-alpha-tubulin captured with CFI Apochromat TIRF 100XC

Photo courtesy of Alexey Khodjakov, Ph.D Research Scientist VI / Professor, Wadsworth


#### NAMC image

Mouse embryos, captured with CFI S Plan Fluor ELWD NAMC 20XC objective

HeLa cells expressing BRET-based calcium indicator protein, Nano-lantern (Ca<sup>2+</sup>). Sample courtesy of Prof. Takeharu Nagai, The Institute of Scientific and Industria Nikon's high-precision CFI60 infinity optics, designed for use with a variety of sophisticated observation methods, are highly regarded by researchers for their superb optical performance and solid reliability.

### Apodized phase contrast

Nikon's unique apodized phase contrast objectives with selective amplitude filters dramatically increase contrast and reduce halo artifacts to provide detailed highdefinition images.

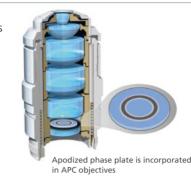

#### **DIC** (Differential Interference Contrast)


Nikon's highly-regarded DIC optics provide uniformly clear and detailed images with high resolution and contrast throughout the magnification range. DIC prisms are individually tailored for each objective lens to provide the highest-quality DIC images for every sample.



DIC prisms matched to individual objectives are mounted in the nosepiece

#### Volume Contrast (Ti2-E)

Volume Contrast imaging utilizes label-free, brightfield images captured at various Z-depths to assemble a phase distribution image. Volume Contrast renders cells easily identifiable as objects for automated counting and area analysis. As this method utilizes brightfield imaging, Volume Contrast enables in-line, non-destructive analysis of cells, suitable for various applications.

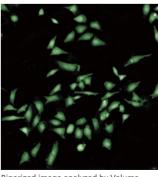


Binarized image analyzed by Volume Contrast

#### External phase contrast (Ti2-E)


The motorized external phase-contrast system enables users to combine phase contrast with epi-fluorescence imaging without compromising fluorescent light transmission by bypassing the need to use phase-contrast objectives. For example, very high NA, liquid immersion objectives can

contrast imaging. Using this external phase contrast system, users can easily combine phase Objective contrast with other imaging modalities, including weakfluorescence imaging such as TIRF and laser tweezer applications.

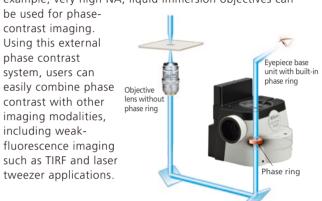


#### NAMC (Nikon Advanced Modulation Contrast)


This is a plastic-compatible, high-contrast imaging technique for unstained, transparent samples such as oocytes. NAMC provides pseudo-three-dimensional images with a shadow-cast appearance. The direction of contrast can be easily adjusted for each sample.

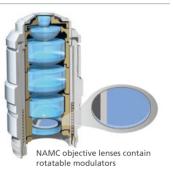


## **Epi-fluorescence**

The Lambda D series objectives, utilizing Nikon's proprietary Nano Crystal Coat technology, are perfect for demanding, low-signal, multi-channel fluorescence imaging that requires high transmission and aberration correction over a wide wavelength range. Combined with new fluorescence filter cubes that offer improved fluorescence detection and stray light countermeasures such as the Noise Terminator, the Lambda D series objectives demonstrate their power in weak signal observations such as single-molecule imaging and even luminescence-based applications.

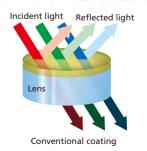

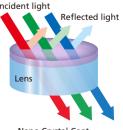

Nano Crystal Coat

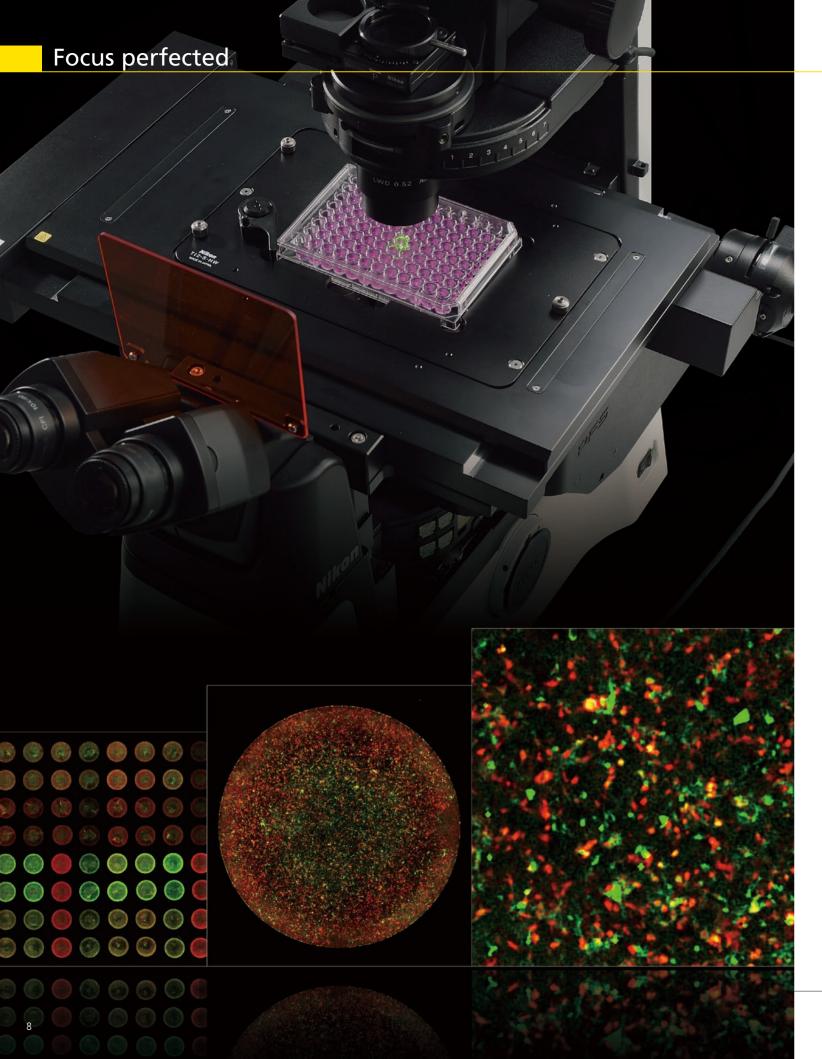

Even the slightest change in temperature and vibrations in the imaging environment can greatly impact focus stability. The Ti2 eliminates focus drift using both static and dynamic measures to enable faithful visualization of the nanoscopic and microscopic world during long time-lapse experiments.

### Mechanically redesigned for ultra-high stability (Ti2-E)

In order to improve the focusing stability, both Z-drive and PFS autofocusing mechanisms have been completely re-designed. The new Z-focusing mechanism is smaller and positioned adjacent to the nosepiece to minimize vibrations. It remains adjacent to the nosepiece even in an expanded (staged-up) configuration, ensuring stability for all applications.



High stability Z-focusing mechanism remains adjacent to the nosepiece even in expanded configurations

The detector portion of the Perfect Focus System (PFS) has been detached from the nosepiece in order to reduce mechanical load on the objective nosepiece. This new design also minimizes heat transfer, which contributes to a more stable imaging environment. Towards this end, the power consumption of the Z-drive motor has also been reduced.


Combined, these mechanical redesigns result in an ultra-stable imaging platform, perfectly suited for single-molecule imaging and super-resolution applications.



### Real time focus correction with the PFS: Simply perfect (Ti2-E)

The Perfect Focus System (PFS) automatically corrects focus drift caused by temperature changes and mechanical vibrations, which can be caused by a variety of factors including the addition of reagents to the sample and multi-position imaging.

The PFS maintains focus by detecting and tracking the position of the coverslip surface in real time. Unique optical offset technology allows users to easily maintain focus at a desired position offset from the cover slip surface. The PFS automatically and continuously maintains focus by means of a built-in linear encoder and high speed feedback mechanism, providing highly reliable images even during long-term, complex imaging tasks

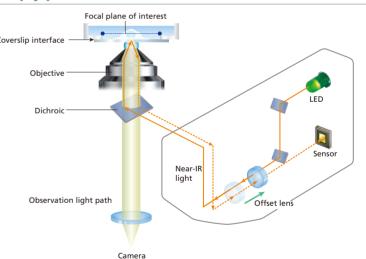

## Water Immersion Dispenser (Ti2-E)

The performance of long-term imaging using the PFS together with water immersion objectives can be increased by using the new Water Immersion Dispenser. The Water Immersion Dispenser automatically applies the appropriate amount of pure water to the tip of an objective, preventing the immersion liquid from drying out and overflowing during experiments. It is compatible with all types of water immersion objectives and helps to stably provide high-resolution, high-contrast and aberration-corrected time-lapse images over long periods of time.



Dual micro nozzle geometry automatically keeps the proper amount of water on the tip of water immersion objectives.



It is no longer necessary to memorize complex microscope alignment and operation procedures. The Ti2 integrates data from sensors to guide you through these steps, eliminating user error and enabling researchers to concentrate on their data.

## Continuous display of microscope status (Ti2-E/A)

A collection of built-in sensors detects and relays status information for a variety of components in the microscope. All of the status information is recorded in the metadata when you acquire images with a computer, so you can easily recall acquisition conditions and/or check for configuration errors.

In addition, a built-in internal camera allows users to view the back aperture, facilitating confirmation of phase ring alignment and extinction cross in DIC. It also provides a laser-safe method for aligning lasers for applications such as TIRF.


Microscope status can be viewed on a tablet and also determined based on status lights on the front of the microscope, enabling status determination in a dark room.





## Guidance for operational procedures (Ti2-E/A)

The Ti2's Assist Guide function provides interactive step-by-step guidance for microscope operation. The Assist Guide can be viewed on a tablet or PC, and integrates real time data from built-in sensors and an internal camera. The Assist Guide is designed to help users through alignment procedures for both experiment setup and troubleshooting.

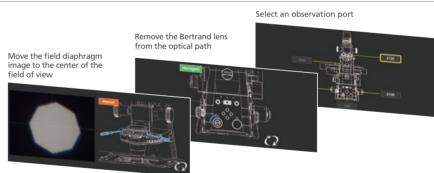

## Automatically detect errors (Ti2-E/A)

The Check Mode allows users to easily confirm, on either a tablet or PC that all the correct microscope components are in place for their chosen observation method. This capability eliminates time and effort normally required for troubleshooting when the desired observation method is not achieved. This functionality is particularly advantageous when multiple users are involved, each with the potential to make unexpected changes to the microscope settings. Custom check procedures can also be pre-programmed.






The Ti2 has been completely redesigned, from the overall body design to the selection and placement of every button and switch, for the ultimate in user experience. The controls are easy to use even in the dark, where the majority of imaging experiments are conducted. The Ti2 provides an intuitive and effortless user interface so researchers can focus on the data and not on microscope control.

### Thoughtfully designed layout for microscope control (Ti2-E/A)

The placement of all of the buttons and switches are based on the type of illumination they control. Buttons that control diascopic observation are positioned on the left side of the microscope and those that control epi-fluorescence observation are on the right side. Buttons that control common operations are on the front panel. This use of zoning provides an easy-to-remember layout, a desirable feature when operating the microscope in a dark room.

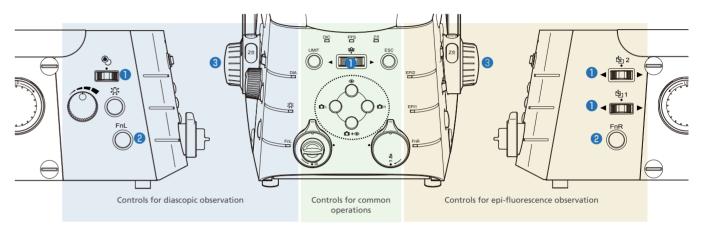

#### 1 Shuttle switch (Ti2-E)

Shuttle switches have been incorporated into the design to control devices such as the epi-fluorescence filter cube turret and objective nosepiece. These types of switches emulate the feel of manually rotating these devices, for intuitive control. Additional functionality can be incorporated into these shuttle switches so that a single switch can operate multiple related devices. For example, the shuttle switch for the epi-fluorescence filter cube turret not only rotates the turret but also opens and closes the fluorescence shutter when the user presses the switch. It is also possible to program these switches to operate a barrier filter wheel and the external phase contrast unit.

#### 2 Programmable Function button (Ti2-E/A)

Conveniently located Function buttons allow customization of the user interface. Users can select from more than 100 functions, including control of motorized devices such as shutters and even signal output to external devices via the I/O port for triggered acquisition. Mode functions, which enable instant changing of observation methods by storing the settings of each motorized device, can also be assigned to these buttons.

#### **3** Focusing knob (Ti2-E)

A focus acceleration button and a PFS engagement button are provided adjacent to the focusing knobs. The two buttons are easily identified by touch because of their different shapes. Focusing speed is automatically adjusted for the objective in use, enabling stress-free operation by maintaining an ideal focusing speed.

## Intuitive control with joystick and tablet (Ti2-E)

The Ti2 joystick not only controls stage movement, but the majority of motorized functions on the microscope, including PFS activity. It can display XYZ coordinates and the status of microscope components, providing an effective means for the user to remotely control the microscope. Motorized functions of the Ti2 can also be controlled from a tablet, connected by wireless LAN to the microscope, providing a versatile graphical interface for microscope control.





System expandability

#### Stage up kit

Allows expansion of the infinity space for incorporation of additional devices such as a second epi-fluorescence filter cube turret, barrier filter wheel, back port unit, and LAPP modules.




#### LAPP system

Various illumination modules, such as TIRF and photostimulation modules, can be flexibly combined to customize your imaging system. Up to five modules can be mounted simultaneously.



### Epi-fluorescence filter cube turret

Compatible with large FOV. A manual intelligent model and high-speed motorized models with motorized shutter are available.

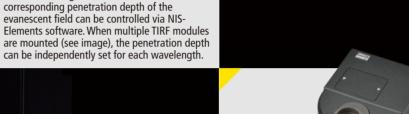

#### Field stop sliders

Two different rectangular-shaped apertures and one round aperture model are available. The rectangular models prevent excitation outside of the imaging area and unintentional photobleaching of samples. They can be removed for ultra-wide FOV imaging.



switch time between adjacent positions) and is compatible with large FOV. It can also be mounted in the infinity space, under the epifluorescence filter cube turret, when the stage is



Control methods are optimized for fast stage

and a piezo stage for Z-stacking are available.

operation. A stage with a built-in encoder

Magnetic specimen holders ensure safe

#### Manual stage


A long travel stroke allows observation of the entire area of a well plate. The travel area can be restricted to ensure safe operation. Various specimen holders are available

Motorized TIRF Module

The incident angle of the laser and

corresponding penetration depth of the

evanescent field can be controlled via NIS-

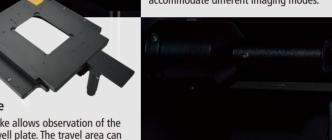

#### Condenser turret

Three types are available: manual, intelligent, and a motorized type that automatically switches between seven positions to accommodate different imaging modes.



The AX high-resolution/AX R ultrafast highresolution confocal system. Multiphoton versions A1 MP+/A1R MP+ are also available.



#### **Auto Correction Collar**

Changes in sample/coverglass thickness, refractive index distribution and temperature can lead to image deterioration. This unique auto correction collar utilizes a harmonic drive and automatic correction algorithm to achieve precise collar adjustment.



#### Fluorescence LED light source

An eco-friendly light source optimized for fluorescence observation. It is a light source that can be directly mounted on the epi-fluorescence module.



#### Oil hydraulic micromanipulator

The MTK-1-N4 achieves simple, easy pipette alignment and smooth pipette movement by means of a four-axis hydraulic joystick. Manufactured by Narishige Lifemed Co., Ltd.



#### Stage top incubator

The STX series precisely maintains temperature at 37.0°C and humidity at more than 95%, and controls CO<sub>2</sub> to enable culturing of cells for more than 1 week.

Manufactured by Tokai Hit Co., Ltd.



#### **Super Resolution Microscopes**

N-SIM S, which achieves live-cell imaging with double the resolution of conventional optical microscopes, and N-STORM, which achieves ten times the resolution of conventional optical microscopes for molecular-level observation, are

# Elements


## A unified acquisition and analysis software platform


Nikon's unified software platform, NIS-Elements, provides acquisition control for basic to advanced imaging systems as well as powerful analysis tools and stunning display capabilities. NIS-Elements can be streamlined for simple, turnkey use and expanded for fully customized, complex experiments such as conditional workflows. NIS-Elements also offers easy-to-use, graphical programming tools such as JOBS and Illumination Sequence for customizing tasks. In addition to Nikon hardware, NIS-Elements controls devices from a vast array of manufacturers to enable the highest level of customization.

## **Device Control**

#### Multidimensional Imaging

Optical Configuration (OC) settings memorize custom observation

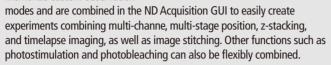

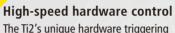

capabilities remove software

callbacks during acquisition routines to maximize imaging speed. The Illumination Sequence module provides a simple, graphical interface for quickly designing and verifying complex, triggered acquisition

#### Graphical programming for custom tasks

For complex or unique image

acquisition needs, the JOBS tool provides an easy-to-use graphical interface for creating custom workflows using drag-and-drop features.

# Control of Third Party Products

NIS-Elements can natively control a vast number of devices from a variety of

manufacturers including high-sensitivity cameras, piezo-devices, light sources, wheels, and National Instruments DAQ devices. This flexibility in hardware control enables custom configurations tailored to individual research needs.


## **Display & Processing**

## **Image Analysis**

#### Automatic measurement

A combination of powerful segmentation tools, morphology

functions, classifiers, and an extensive list of measurement tools for 2D, 3D and timelapse datasets enable users to extract quantitative information from their data with ease. Interactive/manual measurement options are also available.

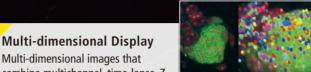

combine multichannel, time-lapse, Z stack, and multi-XY position imaging, as well as image stitching are intuitively displayed in a single window. Powerful volume rendering tools provide unrivalled image quality and interactivity while the

Movie-Maker feature enables users to easily create stunning movies.

#### Real-time measurement

Time measurements can be carried out in real time and visualized

during acquisition. Real-time analysis results can be especially useful for optimizing experiments that rely on drug treatments or ratiometric imaging like FRET or FRAP.




smoothing, and denoising as well as real-time shading correction and image averaging for noise reduction are available. NIS-Elements also offers advanced image arithmetic tools as well as a variety of image projection options such as extended depth of focus (EDF).

### 2D and 3D Object Tracking

NIS-Elements provides powerful tools for identifying and tracking 2D

and 3D objects. Measurements include velocity, acceleration, distance, and direction. A variety of innovative display options are available for communicating and presenting tracking analysis results.

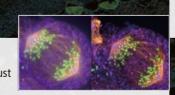

#### Deconvolution Automatic and manual modes, robust

algorithms for noise measurement

Advanced Image Processing Advanced filters for sharpening,

and removal, and enhanced spherical aberration correction are provided to help actualize theoretical resolutions for even confocal images. Both 3D and 2D deconvolution options are available.



### **Custom Analysis Routines**

The General Analysis (GA) module provides an easy method for


creating custom analysis routines that combine image processing and measurement. Routines created in GA can be saved and recalled and even combined with JOBS to create conditional acquisition workflows that rely on real-time analysis results from the integrated GA routine.









## System Diagram

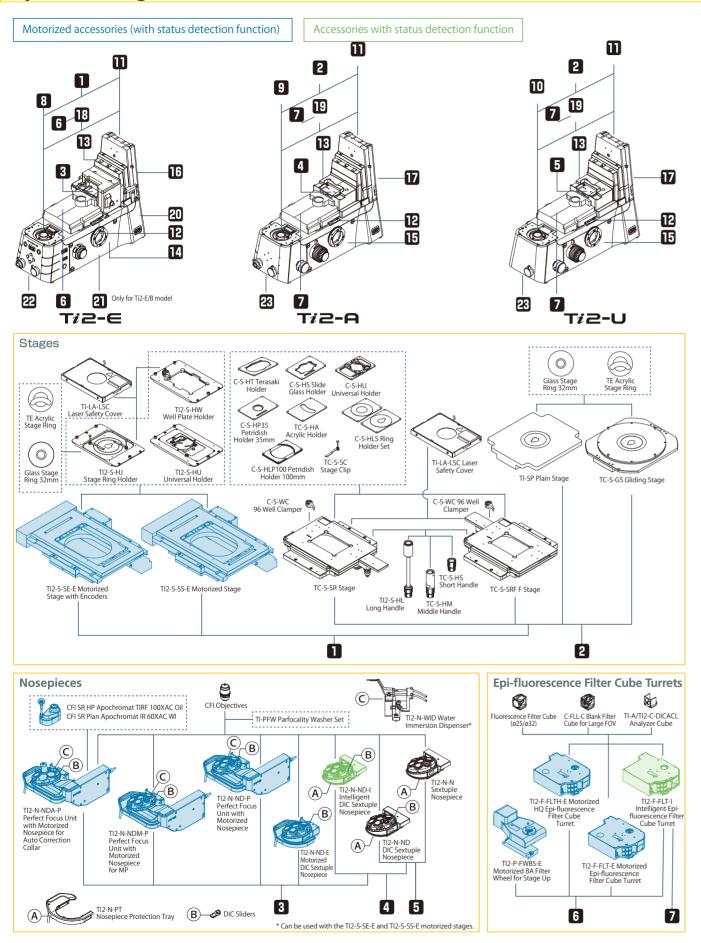


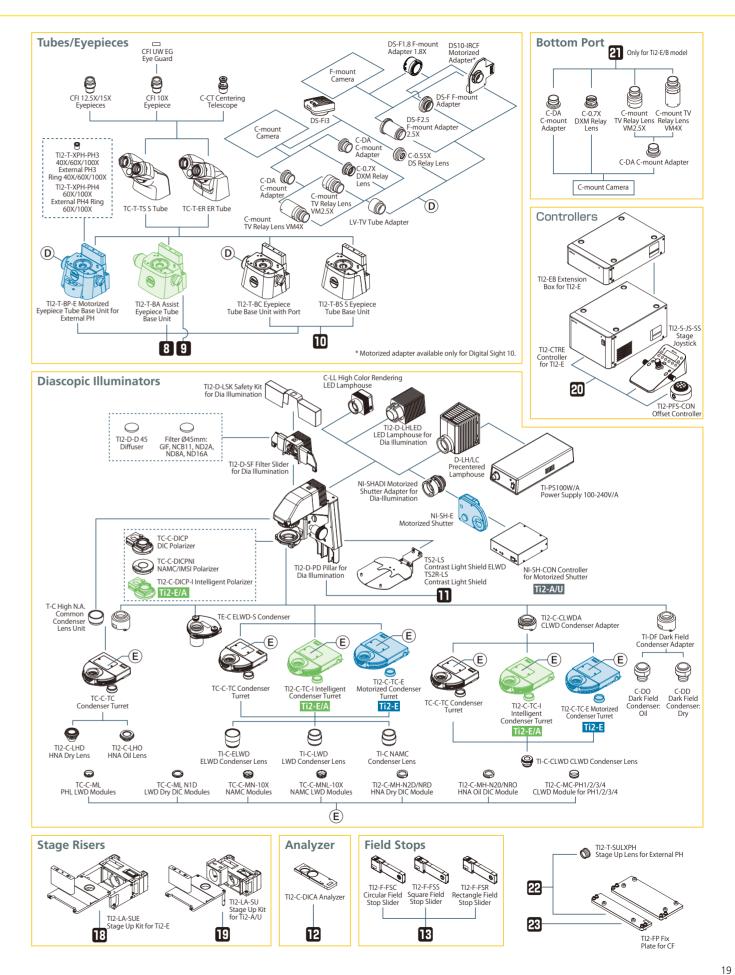

#### Motorized accessories (with status detection function)

Accessories with status detection function

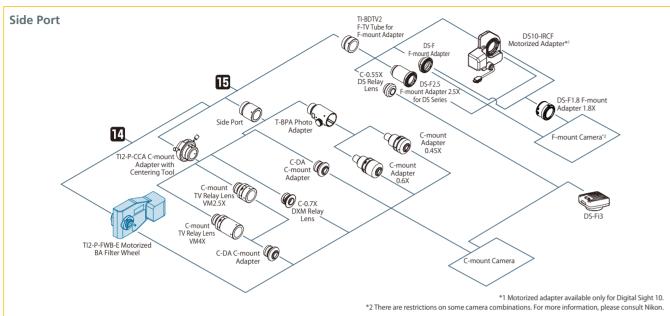

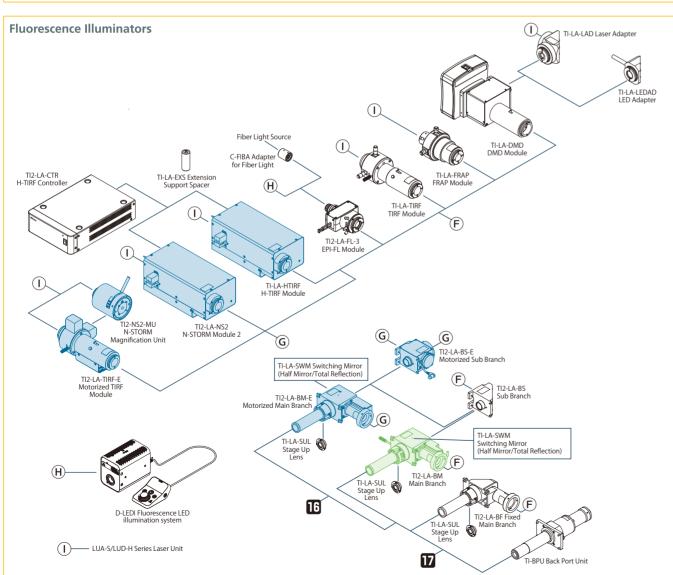

## Objectives

Туре	Model		Immersion	NA	W.D. (mm)	Cover glass	Correction	Spring	DIC	Phase contrast	Fluorescence			PFS
≥			minersion		W.D. (IIIII)	thickness	ring	loaded			UV	Visible light	NIR	113
		4X		0.13	17.20						0	0		
		10X		0.30	16.00	0.17			0		0	0		1
		20X	0.1	0.50	2.10	0.17			0		0	0		_
	2	20XC MI	Oil Glycerin Water	0.75	0.51-0.35 0.51-0.34 0.49-0.33	0-0.17	1	1	0		0	0		
		40X		0.75	0.66	0.17		1	0		0	0		1
		40X Oil	Oil	1.30	0.24	0.17		<b>√</b> w/stopper	0	EXT PH3-40x	0	0		1
_		60XC		0.85	0.40-0.31	0.11-0.23	1	1	0		0	0		
Plan Fluor	CFI Plan Fluor	60XS Oil	Oil	0.50-1.25	0.22	0.17		1	0	EXT PH3-60x	0	0		
Plan		100X Oil	Oil	1.30	0.16	0.17		<b>√</b> w/stopper	0		0	0		1
		100XS Oil	Oil	0.50-1.30	0.16	0.17		✓	0		0	0		
		DL 4XF		0.13	16.50	1.20				PHL	0	0		
		DLL 10X		0.30	16.00	0.17				PH1	0	0		1
		DL 10XF		0.30	15.20	1.20				PH1	0	0		1
		DLL 20X		0.50	2.10	0.17				PH1	0	0		1
		DLL 40X		0.75	0.66	0.17		1		PH2	0	0		1
		DLL 100X Oil	Oil	1.30	0.16	0.17		<b>√</b> w/stopper		PH3	0	0		1
		ADH 100X Oil	Oil	1.30	0.16	0.17		<b>√</b> w/stopper		PH3	0	0		1
		LWD 20XC		0.70	2.30-1.30	0-1.80	1		0		0	0		1
		ELWD 20XC		0.45	8.20-6.90	0-2.00	1		0		0	0		1
		ELWD 40XC		0.60	3.60-2.80	0-2.00	1		0		0	0		1
_		ELWD 60XC		0.70	2.60-1.80	0.10-1.30	1		0		0	0		
Plan Fluo	CELS Blan Elver	LWD ADM 20XC		0.70	2.30-1.30	0-1.80	1			PH2	0	0		1
Plan	CFI S Plan Fluor	ELWD ADM 20XC		0.45	8.20-6.90	0-2.00	1			PH1	0	0		1
S		ELWD ADM 40XC		0.60	3.60-2.80	0-2.00	1			PH2	0	0		1
		ELWD ADL 60XC		0.70	2.60-1.80	0.10-1.30	1			PH2	0	0		
		ELWD NAMC 20XC		0.45	8.20-6.90	0-2.00	1				0	0		
		ELWD NAMC 40XC		0.60	3.60-2.80	0-2.00	1				0	0		
		4X		0.20	15.50	_					◎ 340	0		1
		10X		0.50	1.10	0.17		1	0		◎ 340	0		1
Fluor		20X		0.75	1.00	0.17		1	0		◎ 340	0		1
Super Fluor	CFI Super Fluor	40XC		0.90	0.34-0.26	0.11-0.23	1	1	0		◎ 340	0		
S		40X Oil	Oil	1.30	0.19	0.17		✓w/stopper	0		◎ 340	0		1
		100XS Oil	Oil	0.50-1.30	0.20	0.17		1			◎ 340	0		
	CFI Plan Apochromat	Lambda D 2X		0.10	8.50	0-0.17					© CF	0	0	
		Lambda D 4X		0.20	20.00	0-0.17					0	0	0	1
		Lambda D 10X		0.45	4.00	0.17			0		0	0	0	1
		Lambda D 20X		0.80	0.80	0.17		1	0		0	0	0	1
		Lambda D 40XC		0.95	0.21	0.11-0.23	1	1	0		© CF	0	0	1
		Lambda D 60X Oil	Oil	1.42	0.15	0.17		1	0	EXT PH3-60x	0	0	0	1
		Lambda D 100X Oil	Oil	1.45	0.13	0.17		1	0	EXT PH3-100x	0	0	0	1
		Lambda S 10X		0.45	4.00	0.17			0		•	0	0	1
omat		Lambda S 25XC Sil	Silicone Oil	1.05	0.55	0.11-0.23	1		0		•	0		1
Plan Apochromat		Lambda S 40XC Sil	Silicone Oil	1.25	0.30	0.13-0.21 (23°C)	1		0		•	0		/
an Ap						0.15-0.23 (37°C)				EVE DUO CO.				
E		VC 60XC WI**	Water	1.20	0.31-0.28	0.15-0.18	/	/	0	EXT PH3-60x	0	0		/
		IR 60XC WI**  10XC Glyc	Water Oil Glycerin	0.50	0.18-0.16 2.00***	0.15-0.19	1	/	0	EXT PH3-60x	•	0	0	1
	CELLID DI C	VC 400V 0'	Water	4	0.15	0.77				EVE DIVE 12				
	CFI HP Plan Apochromat	VC 100X Oil	Oil	1.40	0.13	0.17		1	0	EXT PH3-100x	•	0		/
	CFI SR Plan Apochromat	IR 60XC WI**	Water	1.27	0.18-0.16	0.15-0.19	1		0	EXT PH3-60x	•	0	0	/
	CELED LID DI-	IR 60XAC WI*,**	Water	1.27	0.18-0.16	0.15-0.19	1		0	EXT PH3-60x	•	0	0	1
	CFI SR HP Plan Apochromat	Lambda S 100XC Sil	Silicone Oil	1.35	0.31-0.29 (23°C) 0.30-0.28 (37°C)	0.15-0.19	1		0		0	0		1
	CFI Apochromat	LWD Lambda S 20XC WI**	Water	0.95	0.99-0.90	0.11-0.23	1		0			0	0	1
		LWD Lambda S 40XC WI**	Water	1.15	0.61-0.59	0.15-0.19	1		0	EXT PH3-40x	0	0		1
		Lambda S 40XC WI**	Water	1.25	0.20-0.16	0.15-0.19	1	1	0	EXT PH3-40x	0	0		1
Apochromat		TIRF 60XC Oil	Oil	1.49	0.16-0.10 (23°C) 0.13-0.07 (37°C)	0.13-0.19 (23°C) 0.15-0.21 (37°C)	1		0	EXT PH4-60x	•	0		1
Apoc		TIRF 100XC Oil	Oil	1.49	0.16-0.10 (23°C) 0.15-0.09 (37°C)	0.13-0.19 (23°C) 0.14-0.20 (37°C)	1		0	EXT PH4-100x	•	0		1
4	CELSR HP Anochromat	TIDE 100VC Oil	Oil	1.49	0.16-0.10 (23°C)	0.13-0.19 (23°C)	1		0	EXT PH4-100x	•	0		1
1	CFI SR HP Apochromat	TIRF 100XC Oil	Oil	5	0.15-0.09 (37°C) 0.16-0.10 (23°C)	0.14-0.20 (37°C) 0.13-0.19 (23°C)	•							+

<sup>○:</sup>recommended for best results

<sup>•:</sup> possible with visible light that has a longer wavelength than excitation light used for DAPI

CF: confocal imaging is possible at 488 nm and above

EXT: external phase contrast modules 340: high transmittance with an ultraviolet wavelength range of up to 340nm

<sup>\*</sup> Compatible with Auto Correction Collar \*\* Compatible with Auto Correction Collar

\*\* Compatible with Water Immersion Dispenser

\*\*\* 5.50 for upright microscopes

## **Specifications**


		ECLIPSE Ti2-E, Ti2-E/B*1	ECLIPSE Ti2-A	ECLIPSE Ti2-U				
	Optical system	Infinity-corrected CFI60						
	Field number*2	22 with C-mount, 25 with F-mount						
	Intermediate Magnification switching	Manual switching of 1.0x/1.5x (exchangeable from 1.5x to 2.0x)						
	Intermediate Magnification switching	Status detection		_				
	Bertrand lens	Manual in/out, and manual focus, Status detection		_				
Main body	Output port	4 Motorized positions Eyepiece 100%, left 100%, right 100%, eyepiece 20%/left 80% (Ti2-E/B: bottom 100%)	100%, right 100%, option (to 1% or eyepiece 20%/right 80%)					
		Can add ports by use of back port unit and/or choice of tube base	add ports by use of back port unit and/or choice of tube base unit*3					
	Focusing unit	Motorized drive, Coarse/fine focus changeover, 10mm stroke, Minimum increments: 0.01µm, 0.02µm (with encoder control)  Manual drive, Coarse/fine focusing knob, 10mm						
	Stage up	Available*4						
	Binocular tube	Binocular S tube TC-T-TS (field number 22), Ergonomic ER tube TC-	-T-ER (field number 22)					
Tube body	Motorized eyepiece tube base unit for external PH (TI2-T-BP-E)	Camera port (field number 16), Motorized PH turret with 4 motorized positions		_				
	Assist eyepiece tube base unit (TI2-T-BA)	Assist camera (field number 22), Status detection		_				
	Eyepiece tube base unit with port (TI2-T-BC)	Camera port (field number 16)	_	Camera port (field number 16)				
	Pillar for transmitted illumination (TI2-D-PD)	Condenser vertical stroke: 66mm, Backward tilting up to 25 degre 2 filter slot positions (4 filter position option is also available with						
Transmitted illumination	LED Lamphouse for dia illumination (TI2-D-LHLED)	High power LED						
	High Color Rendering LED Lamphouse (C-LL)	High color rendering LED						
	Precentered Lamphouse (D-LH/LC)	100W halogen bulb (pre-centered)						
	Motorized condenser turret (TI2-C-TC-E)	7 motorized positions (Ø37mm x4, Ø39mm x3), LWD/ELWD/CLWD/NAMC condenser lenses are supported		_				
	Intelligent condenser turret (TI2-C-TC-I)	7 manual positions (Ø37mm x4, Ø39mm x3), Status detection, LWD/ELWD/CLWD/NAMC condenser lenses are supported						
Condenser	Condenser turret (TC-C-TC)	7 manual positions (Ø37mm x4, Ø39mm x3), LWD/ELWD/CLWD/NAMC condenser lenses are supported						
	ELWD-S condenser turret (TE-C)	4 manual positions, With ELWD condenser lens (O.D.=65mm, NA=0.3)						
	Condenser lens	LWD (O.D.=30mm, NA=0.52), ELWD (O.D.=75mm, NA=0.3), CLWD (O.D.=13mm, NA=0.72), HNA dry (O.D.=5mm, NA=0.85), HNA oil (O.D.=2mm, NA=1.4), NAMC (O.D.=44mm, NA=0.4)						
	Motorized stage (TI2-S-SE-E, TI2-S-SS-E)	Stroke X: ±57mm, Stroke Y: ±36.5mm, Max drive speed: approx. 25mm/sec, Magnetic sample holder		_				
Stage	Stage (TC-S-SR, TC-S-SRF)	Stroke X: ±57mm, Stroke Y: ±36.5mm, Adjustable stroke range (3 levels) with adjusting pin, Long/middle/short hand options available						
	Gliding stage (TC-S-GS)	Stroke Ø20mm						
	Perfect Focus Unit with motorized nosepiece for Auto Correction Collar (TI2-N-NDA-P)	5 motorized positions, Simple waterproof structure		_				
Nosepiece	Motorized DIC sextuple nosepiece (TI2-N-ND-E) Perfect Focus Unit with motorized nosepiece (TI2-N-ND-P) Perfect Focus Unit with motorized nosepiece for MP (TI2-N-NDM-P)	6 motorized positions, Simple waterproof structure		_				
	Intelligent DIC sextuple nosepiece (TI2-N-ND-I)	6 manual positions, Status detection, Simple waterproof structure		_				
	Sextuple nosepiece (TI2-N-N), DIC sextuple nosepiece (TI2-N-ND)	6 manual positions, Simple waterproof structure						
Epi-fluorescence	Motorized epi-fluorescence filter cube turret (TI2-F-FLTH-E, TI2-F-FLT-E)	6 motorized positions, Motorized shutter		_				
filter cube turret	Intelligent epi-fluorescence filter cube turret (TI2-F-FLT-I)	6 manual positions, Manual shutter, Status detection*5						
Filter wheel/shutter	Motorized BA filter wheel (TI2-P-FWB-E)	7 motorized positions, High speed mode: 50ms, Low vibration mode: 100ms (movement time between adjacent positions)		_				
	Motorized shutter (NI-SH-E)*6	12ms to open/close						
Epi-fluorescence	EPI-FL module (TI2-LA-FL-3)	Supports D-LEDI Fluorescence LED; includes 2-position filter slider and aperture diaphragm In combination with the C-FIBA Adapter for fiber light, fiber illumination can be used						
attachment	Field stop slider	Circular (TI2-F-FSC), rectangular (TI2-F-FSR), square (TI2-F-FSS) a	perture options					
Control unit	Controller, display device	Stage joystick (TI2-S-JS-SS), Tablet	Tablet	_				
Control utill	Controller for TI2-E (TI2-CTRE)	USB/LAN interface, I/O function		_				
Operating environm	nental	Temperature: 0°C+40°C, Humidity: 60% RH max. (at +40°C, no co	ondensation), Indoor use	only				


Motorized accessories have a status detection function

## **Dimensional Diagram**

## T*i* 2-€


Double layer configuration with an Epi-FL module and a FRAP module

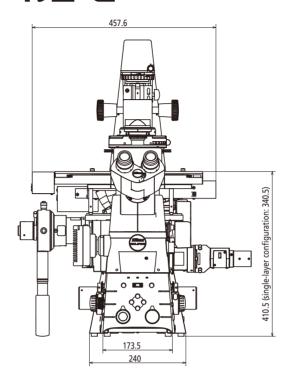

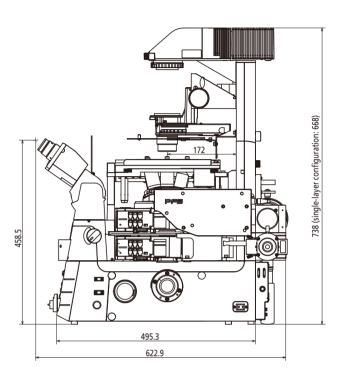

## Ti2-A/U

Single layer configuration with an Epi-FL module

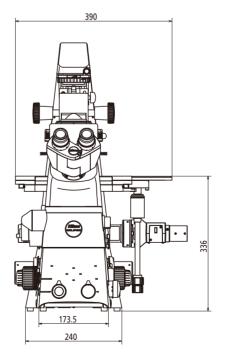

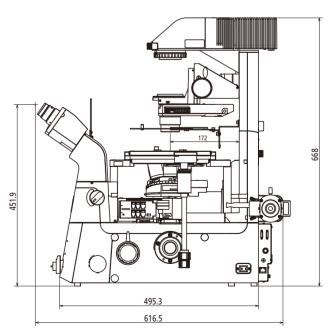


Illustration is of Ti2-A Unit: mm

<sup>\*1</sup> Motorized model with a bottom port

<sup>\*2</sup> Limitations apply based on objective and filter cube choice, stage-up configuration, and illumination module, etc.
\*3 Tube base units with a port cannot be used with Ti2-A

<sup>\*4</sup> Stage up kit is required. Please contact Nikon.
\*5 Status detection cannot be used when attached to the Ti2-U
\*6 NI-SH-CON Controller for Motorized Shutter is required for use with the Ti2-A/Ti2-U

Enter the "Microscopy University" on the web

and discover a whole new world. https://www.microscopyu.com

The Digital Sight series and Imaging Software NIS-Elements are not for clinical diagnostic use.

Specifications and equipment are subject to change without any notice or obligation on the part of the manufacturer. December 2023 © 2018-23 NIKON CORPORATION



TO ENSURE CORRECT USAGE, READ THE CORRESPONDING MANUALS CAREFULLY BEFORE USING YOUR EQUIPMENT.

Monitor images are simulated.

Company names and product names appearing in this brochure are their registered trademarks or trademarks.

N.B. Export of the products' in this brochure is controlled under the Japanese Foreign Exchange and Foreign Trade Law. Appropriate export procedure shall be required in case of export from Japan.

\*Products: Hardware and its technical information (including software)



#### **NIKON CORPORATION**

Shinagawa Intercity Tower C, 2-15-3, Konan, Minato-ku, Tokyo 108-6290, Japan phone: +81-3-6433-3705 fax: +81-3-6433-3785 https://www.healthcare.nikon.com/en/

#### Nikon Instruments Inc.

1300 Walt Whitman Road, Melville, N.Y. 11747-3064, U.S.A. phone: +1-631-547-8500; +1-800-52-NIKON (within the U.S.A. only) . fax: +1-631-547-0299

https://www.microscope.healthcare.nikon.com/

#### Nikon Europe B.V.

Stroombaan 14, 1181 VX Amstelveen, The Netherlands phone: +31-20-7099-000

https://www.microscope.healthcare.nikon.com/en\_EU/

#### Nikon Precision (Shanghai) Co., Ltd.

CHINA phone: +86-21-6841-2050 fax: +86-21-6841-2060 (Beijing branch) phone: +86-10-5831-2028 fax: +86-10-5831-2026 (Guangzhou branch) phone: +86-20-3882-0550 fax: +86-20-3882-0580 https://www.nikon-precision.com.cn/

#### Nikon Canada Inc.

CANADA phone: +1-905-625-9910 fax: +1-905-602-9953

Nikon France, Succursale de Nikon Europe B.V. FRANCE phone: +33-1-4516-4516

Nikon Deutschland, Zweigniederlassung der Nikon Europe B.V.

GERMANY phone: +49-211-9414-888

Nikon Italy, Branch of Nikon Europe B.V. ITALY phone: +39-055-300-9601

Nikon Europe B.V., Amstelveen, Zweigniederlassung Schweiz (Egg/ZH)

SWITZERLAND phone: +41-43-277-2867

Nikon UK, Branch of Nikon Europe B.V. UNITED KINGDOM phone: +44-208-247-1717

Nikon Österreich, Zweigniederlassung der Nikon Europe B.V.

AUSTRIA phone: +43-1-972-6111

Nikon Singapore Pte. Ltd.

SINGAPORE phone: +65-6559-3651 fax: +65-6559-3668

Nikon Australia Pty Ltd AUSTRALIA phone: +61-2-8767-6900

Nikon Instruments Korea Co., Ltd. KOREA phone: +82-2-6288-1900 fax: +82-2-555-4415

NIKON INDIA PVT. LTD.

INDIA phone: +91-124-4688-500

