Inverted Routine Microscope ECLIPSE Ts2



# ECLIPSE Ts2

Inverted Routine Microscope





# Do more than before

# Fits in every Laboratory — Simple to Use & Compact

### Easy to work with

#### Efficient and comfortable observation

LED light sources and built-in Diascopic/Epi-fluorescence illumination systems reduce setup time and allow users to concentrate on their research

#### New streamlined operation

Control buttons on the Ts2 microscope are intuitively located for a streamlined workflow. Commonly used controls such as the on/off and diascopic/epi-fluorescence switching buttons are located at the front panel for easy reach. Buttons pertaining to either diascopic or epi-fluorescence control are zoned to the left and right sides of the microscope body, respectively, to eliminate confusion and improve workflow efficiency.



#### Faster, brighter images with LED illumination

LED light source is alignmentfree resulting in faster setup and consistent results. LEDs also eliminate frequent bulb replacements, saving the user time and money. Moreover, The new Contrast Shield (optional) provides high signal-to-noise fluorescence observation even in brightly lit culture rooms.



#### Easy-to-use Mechanical Stage

The high performance mechanical stage (optional) can accommodate a wide range of flasks and cell culture chambers. The new stage design also directly accommodates micro plates. In addition, the sample holder is easily removed to accommodate large flasks.




## Compact, streamlined body for efficient observation

#### LED-based high-quality diascopic and epi-fluorescence observation

Two models are available to meet your needs: a diascopic illumination model, the Ts2, and an epifluorescence illumination model, the Ts2-FL. High-intensity LED sources are employed for both diascopic

and epi-fluorescence illumination. The built-in fly-eye lens ensures uniform brightness across the entire field of view. A wide range of wavelengths is available to choose from for Epi-fluorescence LED illumination.

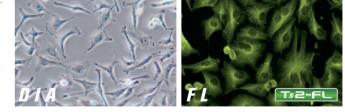

#### Compact and highly stable body


Illumination modules including the epi-fluorescence light source have been seamlessly incorporated into the microscope main body, resulting in a compact and simple design form that's also durable. The compact structure is also vibration-resistant to provide highly stable sample observations

The camera port is located on the side of the microscope to provide unimpeded viewing of the stage even when placed inside a culture hood.







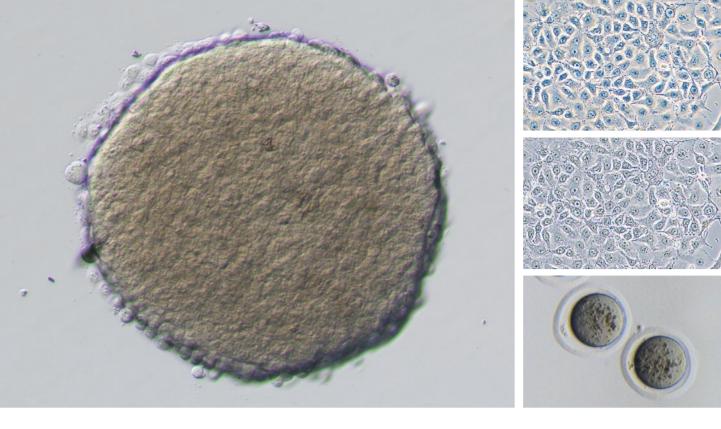

# Do more than before — DIA




# Highly optical performance with diascopic observation


#### Diascopic observation with high-intensity LED (Eco-illumination)

Eco-illumination provides high-intensity light suitable for phase contrast observation. With the built-in fly-eye lens, uniform brightness is provided across the entire field of view. LEDs are an environmentally friendly, low-power-consumption light source. Eco-illumination provides a long lifetime of 60,000 hours and reduces the frequency of lamp replacement.

#### Phase contrast observation

Phase contrast is an optical contrasting technique that typically utilizes a phase contrast objective lens and condenser annulus. The use of a highintensity LED light source results in clear images even at high magnifications

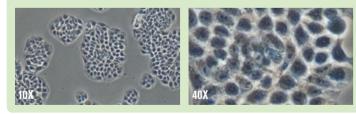

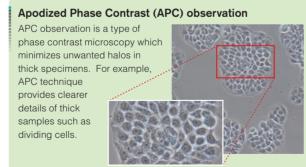


#### New contrasting technique, "Emboss Contrast"


Emboss Contrast is a cost-effective optical technique which does not require costly optics. Utilizing just a bright-field objective lens and two contrast sliders, Emboss Contrast provides pseudo-three dimensional and glare-free images for thick specimens such as iPS cells which would normally suffer from halos with conventional phase contrast methods. Additionally, Emboss Contrast is compatible with both glass and plastic culture chambers, making it a very versatile observation technique.


#### **Comparison of new Emboss Contrast** and Nikon Advanced Modulation Contrast



NAMC / 20X

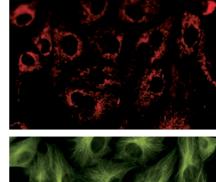

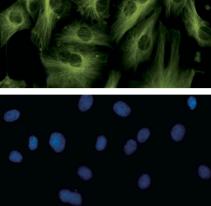

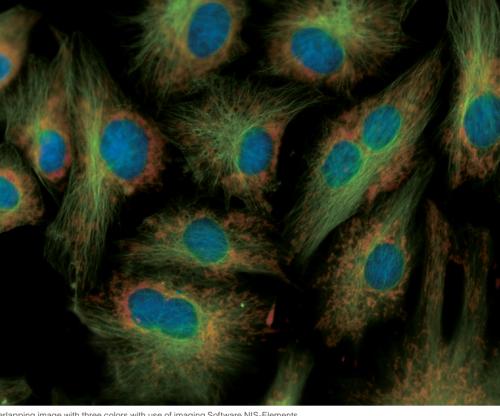

Overlapping image with three colors with use of imaging Software NIS-Elements

# Do more than before — FL Epi-fluorescence observation made easy with LED

Fly-eye lens for uniform illumination

With a built-in fly-eye lens, uniform brightness is provided across the entire field of view.

#### Accurately reproduce illumination power every time


The Ts2 can be configured with up to three fluorescent filter cubes. The illumination power previously defined by the user is replicated when the same wavelength is used again, thus eliminating the need for manual adjustment of light intensity when switching between wavelengths. The Ts2 also incorporates a noise terminator mechanism which allows high signal-to-noise fluorescent images to be captured.





#### High S/N epi-florescence observation in bright rooms

The new Contrast Shield accessory (optional) blocks room light, providing an easy and cost-effective method for achieving high signalto-noise fluorescence observation in a brightly lit culture room.

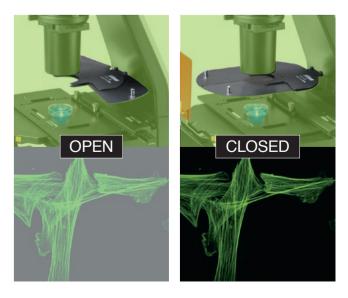

#### Accessories

#### Camera Port

Optional camera port is available for image capturing. Digital Cameras utilizing C- or F-mounts can be attached.



#### Emboss Contrast Slider

Both condenser-side slider and eyepiece-tube-side slider are available. Contrast modules for 10X, 20X, 40X objective lenses are arranged on the same slider so switching between magnifications is easily achieved by simply sliding the contrast slider.



#### ThermoPlate<sup>®</sup> TPi-TS2X (for the Mechanical Stage)

ThermoPlate<sup>®</sup> TPi-TS2X provides accurate and stable temperature control for the specimen from room temperature to 50 degrees Celsius. Proprietary treatment methods ensure that the glass surface of the Termo Plate is breakage-free.

Manufacturer: TOKAI HIT Co.,Ltd.



### Cameras for microscopes


All cameras of the digital sight series can be directly connected to a PC via a fast USB3.0 interface.

\*The optional camera port is required to attach the digital camera to the microscope.





C-mount CMOS Camera				F-mount CMOS Camera
Microscope camera Digital Sight 1000 2.0 megapixel Color Full HD	Nikon Papar spin Mos	Microscope camera DS-Fi3 5.9 megapixel Color High-resolution		Microscope camera Digital Sight 10 23.9 megapixel Color/Monochrome High-resolution
Equipped with a 2 megapixel CMOS image sensor, the Digital Sight 1000 can display, capture and save full HD, 1920x1080 pixel images at 30 fps without using a PC.		A high-definition 5.9-megapixel color CMOS image sensor captures fine-textured images in faithful color. For image acquisition, NIS- Elements imaging software is required.		Achieves color / monochrome switching shooting with a single camera. You can quickly shoot 6K high-definition images in one shot.
Frame rate	30 fps (1920×1080	0) 15 fps (2880 × 2048), 30 fps (1440 × 1024)		9 fps (6000 × 3984), 66 fps (1920 × 1080)
Max recordable pixels	1920×1080	2880 × 2048		6000 × 3984
Imaging sof NIS Advanced Solutions for your lim for a tablet PC	ftware control of cameras,	talling NIS-Elements L on a tablet PC enables s Digital Sight 1000/DS-Fi3/Digital Sight 10 micro live image display, and image acquisition. ation about compatible tablet PCs, contact Nikon.	<u> </u>	

# System diagram

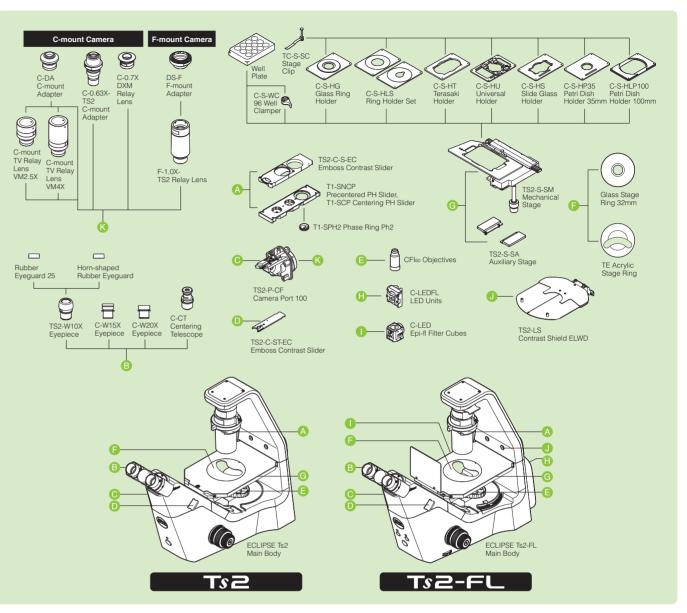

## Dimensions (Unit: mm)

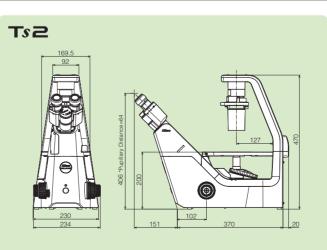

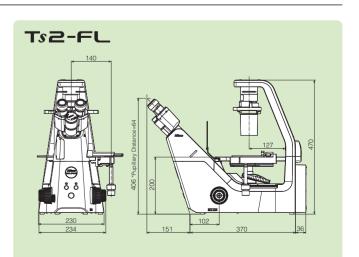

#### T=P / T=P-FL Specifications

192 / 192-FL Speci	ICALIONS			
	Тя2	Ts2-FL		
Optical System	CFIeo Infinity Optical System			
Observation method	Brightfield, Apodized Phase Contrast <sup>*1</sup> , Phase Contrast, Emboss Contrast <sup>*2</sup>	Brightfield, Apodized Phase Contrast*1, Phase Contrast, Emboss Contrast*2, Epi-Fluorescence		
Illumination Diascopic illumination	High luminescent white LED illuminator (Eco-illumination), Built-in Fly eye lens			
Episcopic illumination	—	LED illuminator, built-in Fly eye lens, Can be configured with up to 3 different fluorescence LED units; available wavelengths: 385, 455, 470, 505, 525, 560, 590, 625 nm		
Tube	Inclination: 45 degree, Pupillary distance: 50 - 75 mm, Siedentopf type, Attachable camera port, Eyepiece/Port: 100/0:0/100			
Eyepiece (F.O.V.)	10X (22), 15X (16), 20X (12.5)			
Focusing	Via nosepiece up/down movement, Stroke (manual): Up 7 mm down 1.5 mm Coarse stroke: 37.7 mm per rotation, Fine stroke: 0.2 mm per rotation, Coarse motion torque adjustable			
Nosepiece	Quintuple nosepiece			
Condenser	ELWD Condenser (NA 0.3, W.D. 75 mm)			
Slider	<ul> <li>Precentered or Centering PH Slider, 10X, 20X, 40X Objectives available for phase contrast</li> <li>Emboss Contrast sliders (both condenser-side slider and eyepiece-tube-side slider must be mounted), 10X, 20X, 40X objectives available for Emboss Contrast</li> </ul>			
Stage	<ul> <li>Plain Stage, stage size: 170(X)×247(Y) mm, With 2 Acrylic Types of Stage Ring</li> <li>Mechanical stage (optional), stroke:126(X)×78(Y) mm, Accepts 5 types of micro-testplate, well clamper and stage clip</li> </ul>			
Holder	C-S-HP35 Petridish Holder 35 mm     C-S-HLP100 Petridish Holder 100 mm     C-S-HT Terasaki Holder for Terasaki holder and ø65 dish     C-S-HS Slide Glass Holder for glass slides, ø54 dish and hemocytometer     C-S-HU Universal Holder for Terasaki plate holder, glass slide, ø35-65 dish and hemocytometer     C-S-HG Glass Ring Holder     C-S-HLS Ring Holder Set			
Epi Fluorescence attachment	-	Epi-fluorescence filter turret (with main body), Filter cubes with noise terminator mechanism Configure with up to 3 Epi-fluorescence filter cubes, Additional positions for bright-field observation, Attachable Contrast Shield (optional)		
Dimensions	236(W)×548(D)×471(H) mm	236(W)×564(D)×471(H) mm		
Weight (approx.)	13kg	14.5kg		
Rated Voltage/Electric Current	100 V-240 VAC±10 %, 50/60 Hz, 0.35 A			
Power Consumption	15 W			


\*1 APC (Apodized Phase Contrast) is a type of phase contrast observation with reduced halo, thanks to Nikon's unique lens coating.

\*2 Emboss contrast is Nikon's unique contrast observation method. It provides pseudo-three-dimensional images using focal illumination, which gives high contrast to samples

#### **Related Products** ECLIPSE TS2R/TS2R-FL

A compact inverted microscope for your basic research needs.

Ts2R/Ts2R-FL provides a wide range of observation methods and applications in a compact body that can easily fit in limited laboratory spaces while providing streamlined operation

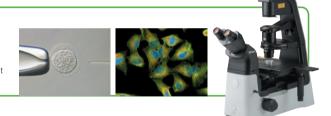

Specifications and equipment are subject to change without any notice or obligation on the part of the manufacturer. July 2024 ©2016-2024 NIKON CORPORATION N.B. Export of the products\* in this brochure is controlled under the Japanese Foreign Exchange and Foreign Trade Law. Appropriate export procedures shall be required in case of export from Japan. \*Products: Hardware and its technical information (including software)

WARNING TO ENSURE CORRECT USAGE, READ THE CORRESPONDING MANUALS CAREFULLY BEFORE USING THE EQUIPMENT.



#### NIKON CORPORATION

Head office 1-5-20, Nishioi, Shinagawa-ku, Tokyo, 140-8601, Japan https://www.healthcare.nikon.com/en/

Manufacturer 471, Nagaodai-cho, Sakae-ku, Yokohama, Kanagawa 244-8533, Japan

ISO 14001 Certified for NIKON CORPORATION

Nikon Instruments Inc. 1300 Walt Whitman Road, Melville, N.Y. 11747-3064, U.S.A. phone: +1-631-547-8500; +1-800-52-NIKON (within the U.S.A. only) fax: +1-631-547-0299

https://www.microscope.healthcare.nikon.com/ Nikon Europe B.V.

Stroombaan 14, 1181 VX Amstelveen, The Netherlands phone: +31-20-7099-000

https://www.microscope.healthcare.nikon.com/en\_EU/

Nikon Precision (Shanghai) Co., Ltd. CHINA phone: +86-21-6841-2050 fax: +86-21-6841-2060 (Beijing branch) phone: +86-10-5831-2028 fax: +86-10-5831-2026 (Guangzhou branch) phone: +86-2-3882-0551 fax: +86-2-3882-0580 https://www.nikon-precision.com.cn/

Nikon Canada Inc.

CANADA phone: +1-905-625-9910 fax: +1-905-602-9953 Nikon France, Succursale de Nikon Europe B.V. FRANCE phone: +33-1-4516-4516

Nikon Deutschland, Zweigniederlassung der Nikon Europe B.V.

GERMANY phone: +49-211-9414-888 Nikon Italy, Branch of Nikon Europe B.V.

ITALY phone: +39-055-300-9601 Nikon Europe B.V., Amstelveen,

Zweigniederlassung Schweiz (Egg/ZH) SWITZERLAND phone: +41-43-277-2867 NIKON UK, Branch of Nikon Europe B.V.

UNITED KINGDOM phone: +44-208-247-171 Nikon Österreich, Zweigniederlassung der Nikon Europe B.V. AUSTRIA phone: +43-1-972-6111

Nikon Singapore Pte Ltd SINGAPORE phone: +65-6559-3651 fax: +65-6559-3668 Nikon Australia Pty Ltd AUSTRALIA phone: +61-2-8767-6900 Nikon Instruments Korea Co., Ltd. KOREA phone: +82-2-6288-1900 fax: +82-2-555-4415 NIKON INDIA PVT. LTD. AUSTRALIA phone: +91-124-4688-500

Code No. 2CE-MPJH-9 (2407) Am